



## NANOTOUCH PRO

CLEANING AND  
DISINFECTION  
SOLUTION

INSTRUCTIONS FOR USE

NON-FLEETING  
PERSISTENT  
DISINFECTION

We recommend that you read the user manual before using this product, keep it during the use, and use the product in accordance with its intended use so that you can get the best efficiency throughout its life.

- Read the user manual before using the product.
- Pay attention to the safety warnings.
- Store the user manual in an easy-to-reach place for later use.
- Read all other documents given with the product.

The below symbol(s) are present in this instruction for use:



Life and property threatening situation

## WHAT IS NANOTOUCH PRO CLEANING AND DISINFECTION SOLUTION?

Nanotouch Pro is a next generation disinfectant that used on all surfaces of the infrastructure such as floors, walls, ceilings, doors, metal and plastic seats, sinks in IVF laboratories; that prevents the development and contamination of fungi, spores, yeast, bacteria, mycoplasma and viruses (including HIV and Hepatitis B) and makes disinfection permanent by binding to the surfaces it is applied to.

### **Nanotouch Pro;**

- ✓ Non-toxic and non-allergic, environmentally friendly,
- ✓ Compatible with many surfaces including aluminum, chrome, nickel, steel, polycarbonate, etc.,
- ✓ Odorless,
- ✓ Alcohol free,
- ✓ VOC free,
- ✓ Does not contain chlorine, aldehyde or phenol, stored at room temperature,
- ✓ Non-corrosive, next-generation disinfectant.

### **USAGE AREAS**

Nanotouch Pro can be safely used in IVF laboratory work surfaces which require hygienic conditions such as ceramic, wood, stone, glass, aluminum, chrome, nickel, steel and polycarbonate or silicone, plastic and polyurethane materials.

### **RECOMMENDED USAGE INTERVAL**

Suitable for daily use.

### **INSTRUCTIONS FOR USE**

#### **Mopping**

1. Take of water into the mopping bucket.
2. Add at least one bottle of Nanotouch Pro (200 ml or optionally increase) to the water.
3. Mix it well.
4. Clean the visible dirt on the floor.
5. Soak the mop in the liquid for at least 1-3 minutes and make sure the liquid is thoroughly absorbed.
6. Spread the liquid in the bucket over the entire floor by mopping.
7. Let it dry.

## Spraying

1. The Nanotouch Pro is optionally diluted up to 1/5 in the spraying device.  
(Add 200 ml to 800 ml of water)
2. Clean the visible dirt on the floor and surface, if there is any.
3. Spray it on all surfaces-to-be-applied from a distance of 45cm without any gaps.
4. Let it dry.



### ATTENTION!

Nanotouch Pro should not be used in any way other than the procedure outlined in this user manual. Do not ingest the product.

### ACTIVE COMPONENT

### SHELF LIFE

### PACKAGING FORMS

Ready-to-use QAC-Si

Minimum 2 years in RT.  
3 months if the package is opened.

200ml, 1L, 5L